

# Analysis of 18 Polycyclic Aromatic Hydrocarbons in Soil Using the QuEChERS Method

Anila I Khan, Thermo Fisher Scientific, Runcorn, Cheshire, UK

## Key Words

PAH, benzofluoranthenes b, j and k, soil, QuEChERS, TraceGOLD TG-17SiIMS

## Abstract

The use of QuEChERS dispersive SPE as a simple, fast, and quantitative sample preparation method is demonstrated for the GC-MS analysis of 18 polycyclic aromatic hydrocarbons (PAHs) in soil. The suitability of the Thermo Scientific™ TraceGOLD™ TG-17SiIMS GC column for the separation of 18 PAHs is also shown. The average recoveries for the spiked 18 PAHs in soil at 1 mg/kg were between 85.0% and 106.7% with relative standard deviations between 0.3% and 2.8% using the original QuEChERS methodology.

## Introduction

Polycyclic aromatic hydrocarbons (PAHs) are some of the most widespread organic pollutants. In addition to occurring naturally in oil, coal, and tar deposits, PAHs are produced by incomplete combustion of fossil fuels and other organic matter. The compounds are of concern because many have been identified as carcinogenic, mutagenic, or teratogenic. Due to these health risks, regulatory agencies, such as the Environmental Protection Agency (EPA), have defined maximum allowable levels of PAHs in the environment. Sensitive analytical methods are essential in the determination of the presence and levels of PAHs.

QuEChERS (Quick, Easy, Cheap, Effective, Rugged and Safe) is a dispersive solid-phase extraction (SPE) technique initially developed for extracting multi-residue pesticides from fruits and vegetables[1]. The technique has now been extended outside of the food safety field to applications such as removing soil matrix from non-pesticide analytes such as PAHs. The advantages of this methodology are speed, ease of execution, minimal solvent requirement, and cost.

The method is:

- Quick – It has high sample throughput. Typically eight samples can be prepared in under 30 min.
- Easy – It requires less handling of extracts than other techniques; i.e., fewer steps are required.



- Cheap – Less sorbent material and less time is needed to process samples compared to other techniques.
- Effective – The simple technique gives high and accurate recovery levels for a range of different compound types.
- Rugged – The method can detect a large number of compounds including non-polar and polar compounds.
- Safe – Unlike other techniques, it does not require the use of any chlorinated solvents. Extraction is typically carried out using acetonitrile, which is both GC and LC compatible.

A sample preparation approach, described in the original QuEChERS procedure [1], was used for extracting 18 PAHs from soil. This is a two-stage process: sample extraction, followed by dispersive SPE.

In the sample extraction stage, the soil sample is sieved to give uniform surface area so that optimal extraction efficiency can be achieved. The soil sample must be at least 80% hydrated for extraction to work; therefore, water must be added to the soil in the extraction tube. This is

then followed by the addition of acetonitrile and a salt mixture composed of magnesium sulfate and sodium chloride. The salt mixture initiates a phase separation between water and organic solvent and the analytes of interest are extracted into the organic phase. The tube is then capped, shaken vigorously, and centrifuged.

The second stage of the QuEChERS method uses dispersive SPE, which involves transferring a portion of the acetonitrile extract to a clean-up tube containing a combination of sorbents for removal of unwanted sample components. The sorbent combination of magnesium sulfate and PSA (primary & secondary amines) silica in the sample clean-up tube removes interfering components from the matrix, thus reducing matrix effects and improving method robustness.

The extraction of 18 PAHs from soil was demonstrated using the original QuEChERS methodology. Six extractions at 1 mg/kg of 18 PAHs spiked into soil were used for recovery experiments. Separation of 18 PAHs was achieved on a Thermo Scientific TraceGOLD TG-17SiIMS GC column. This low-bleed GC column allows for the quantification of critical pairs such as phenanthrene and anthracene. Similarly, it also provides baseline resolution for isobaric PAHs such as benzo[fluoranthene b, j and k.

## Experimental Details

| Consumables                    | Part Number  |
|--------------------------------|--|
| Column:                        | TraceGOLD TG-17SiIMS, 30 m x 0.25 mm x 0.25 µm 26072-1420  |
| Septum:                        | BTO, 17 mm 31303211  |
| Liner:                         | Thermo Scientific™ Splitless FocusLiner™ for 50 mm needle, 5 x 8 x 105 mm 453T2999   |
| Column Ferrules:               | 100% graphite ferrules for Thermo Scientific™ TRACE™ injector 0.1–0.25 mm ID 29053488  |
| Colum Ferrules:                | Graphite/Vespel® for transfer line 0.1–0.25 mm ID 29033496   |
| Injection Syringe:             | 10 µL fixed needle syringe for Thermo Scientific™ TriPlus RSH™ Autosampler 365D0291  |
| Vials and Closures:            | Thermo Scientific 9 mm Wide Opening Screw Thread Vial Convenience Kit, 2 mL Clear Vial with Patch, Blue Polypropylene Closure with Clear PTFE/Blue Silicone Septa 60180-599  |
| Reagents:                      | Fisher Scientific™ LC-MS grade acetonitrile A/0638/17<br>Fisher Scientific LC-MS grade water W/0112/17   |
| QuEChERS Extraction Stage:     | Thermo Scientific™ HyperSep™ Dispersive SPE multipacks, 4000 mg magnesium sulfate, and 1000 mg sodium chloride. Each pack contains 50 metalized pouches with 50 empty centrifuge tubes with plug seal caps 60105-332 |
| QuEChERS Dispersive SPE Stage: | Thermo Scientific HyperSep Dispersive SPE clean-up product, 2 mL tube containing 150 mg magnesium sulfate, 50 mg PSA 60105-203   |

## Preparation of Calibration Standards

The calibration curve was constructed with a mixture of 18 PAHs listed in Table 1 at the following concentrations: 50, 100, 200, 500, 1000, and 2000 ng/mL. To each 1 mL of calibration standard solution, 10 µL of 50 µg/mL mixture of internal standards (IS) containing naphthalene-d<sub>8</sub>, acenaphthene-d<sub>10</sub>, phenanthrene-d<sub>10</sub>, chrysene-d<sub>12</sub>, and perylene-d<sub>12</sub> was added.

| Internal Standard                         | Compound   |
|---|--|
| <b>Napthalene-d<sub>8</sub> (IS 1)</b>    | Napthalene   |
| <b>Acenaphthene-d<sub>10</sub> (IS 2)</b> | Acenaphthylene, Acenaphthene, Fluorene   |
| <b>Phenanthrene-d<sub>10</sub> (IS 3)</b> | Phenanthrene, Anthracene, Fluoranthene, Pyrene, Benzo[a]anthracene   |
| <b>Chrysene-d<sub>12</sub> (IS 4)</b>     | Chrysene, Benzo[b]fluoranthene, Benzo[k]fluoranthene, Benzo[j]fluoranthene, Benzo[a]pyrene, Benzo[e]pyrene |
| <b>Perylene-d<sub>12</sub> (IS 5)</b>     | Indeno[1,2,3-cd]pyrene, Dibenzo[a,h]anthracene, Benzo[g,h,i]perylene                                       |

Table 1: PAHs quantified against the internal standards used for calibration and recovery experiments

## Sample Preparation

Samples were prepared as described in Figure 1.

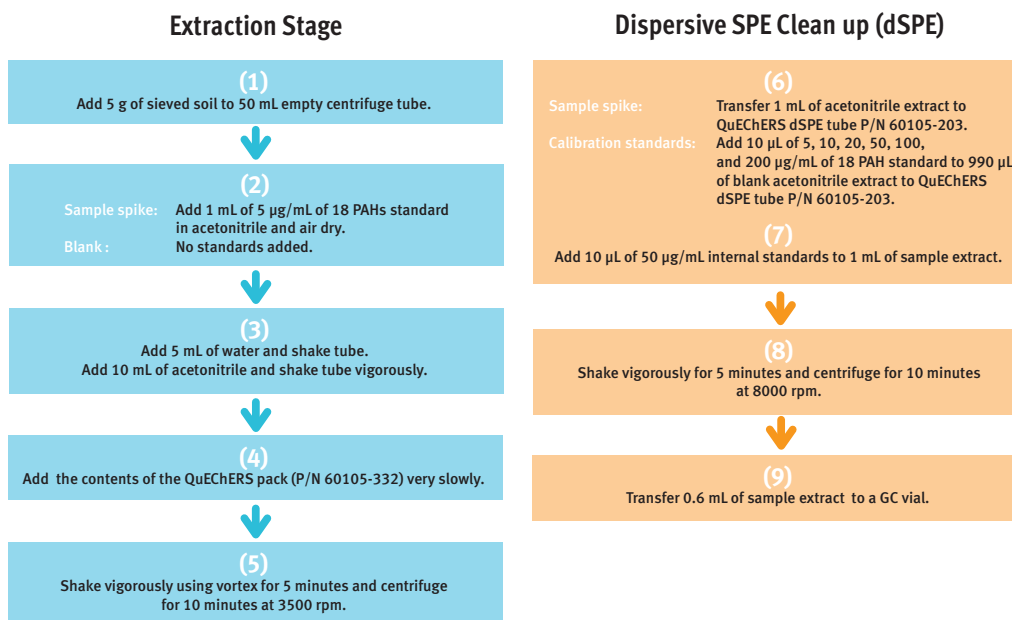


Figure 1: Sample preparation illustrating QuEChERS protocol

## Separation Conditions

|                            |  |
|----------------------------|--|
| Instrumentation:           | Thermo Scientific TRACE GC Ultra                             |
| Carrier Gas:               | Helium   |
| Column Flow:               | 1.2 mL/min, constant flow                                    |
| Oven Temperature:          | 90 °C (1.0 min), 30 °C/min, 250 °C, 4 °C/min, 330 °C (5 min) |
| Injector Type:             | Split/Splitless  |
| Injector Mode:             | Split 25:1, 30 mL/min split flow                             |
| Injector Temperature:      | 250 °C   |
| Detector Type:             | Thermo Scientific™ ISQ™ mass spectrometer                    |
| Transfer Line Temperature: | 300 °C   |
| Source Temperature:        | 250 °C   |
| Ionization Conditions:     | EI   |
| Electron Energy:           | 70 eV  |
| SIM Scan Parameters:       | Table 2  |

| Scan window start time (min) | Compound  | <i>m/z</i> Quan ion | Dwell time (sec) |
|------------------------------|---|---------------------|------------------|
| 3.00                         | Napthalene, Napthalene-d <sub>8</sub>                                 | 128, 136            | 0.12             |
| 4.80                         | Acenaphthylene, Acenaphthene, Acenaphthene-d <sub>10</sub>            | 152, 154, 164       | 0.10             |
| 5.90                         | Fluorene  | 166                 | 0.25             |
| 6.60                         | Phenanthrene, Anthracene, Phenanthrene-d <sub>10</sub>                | 178, 188            | 0.12             |
| 8.10                         | Fluoranthene, Pyrene  | 202                 | 0.25             |
| 11.40                        | Benz[a]anthracene, Chrysene, Chrysene-d <sub>12</sub>                 | 228, 240            | 0.12             |
| 15.50                        | Benzo[b]-, Benzo[j]-, Benzo[k]fluoranthene, Benzo[a]-, Benzo[e]pyrene | 252                 | 0.25             |
| 19.55                        | Perylene-d <sub>12</sub>  | 264                 | 0.25             |
| 22.00                        | Indeno[1,2,3-cd]pyrene, Dibenz[a,h]anthracene                         | 276, 278            | 0.12             |
| 24.80                        | Benzo[g,h,i]perylene  | 276                 | 0.25             |

Table 2: SIM scan parameters

## Injection Conditions

|                                     |   |
|-------------------------------------|---|
| Instrumentation:                    | Thermo Scientific TriPlus RSH Autosampler |
| Injection Volume:                   | 1 $\mu$ L                                 |
| Pre- and Post-Injection Dwell Time: | 1.0 s                                     |

## Data Processing

|           |                              |
|-----------|------------------------------|
| Software: | Thermo Scientific™ XCalibur™ |
|-----------|------------------------------|

## Results

The separation of the 18 PAHs was obtained using a low bleed TraceGOLD TG-17SilMS GC column. The analysis was performed in SIM mode. Figure 2 shows the SIM chromatogram of spiked PAHs in soil matrix at 1 mg/kg level.

To assess the method linearity, a calibration curve was constructed for each of the PAHs using the appropriate internal standard (Table 1), The coefficients of determination ( $R^2$ ) between area ratio of sample and internal standard for 18 PAHs were  $>0.997$  (Table 3), demonstrating excellent method linearity.

Six extractions of spiked samples at 1 mg/kg were carried out and recoveries and reproducibility (RSD) measured. Table 3 shows the recoveries and RSDs for the spiked 18 PAHs in soil, which were between 85.0% and 106.7%, with RSDs between 0.3% and 2.8%.

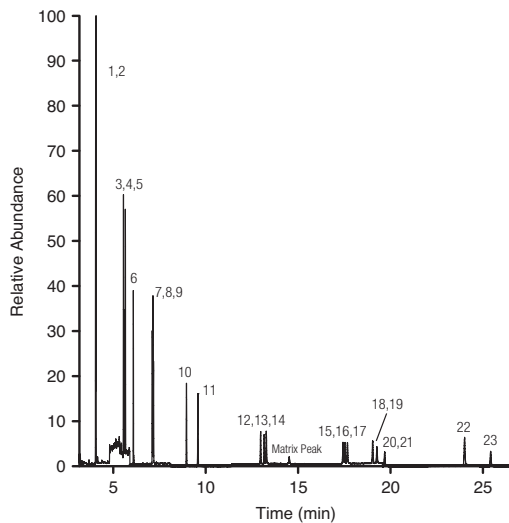


Figure 2a: SIM chromatogram of 18 PAHs in soil matrix solution at 500 ng/mL

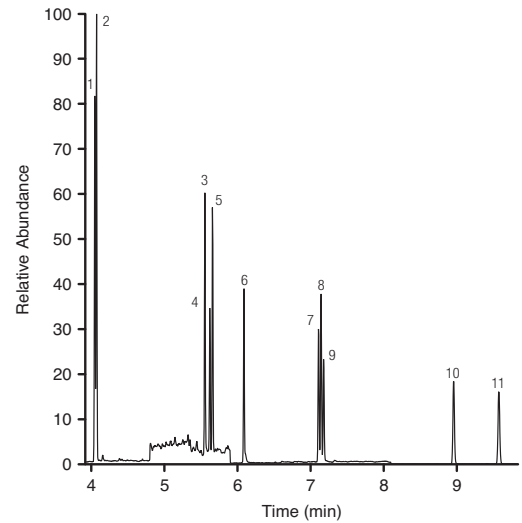


Figure 2b: Expanded SIM chromatogram of 18 PAHs in soil matrix solution at 500 ng/mL

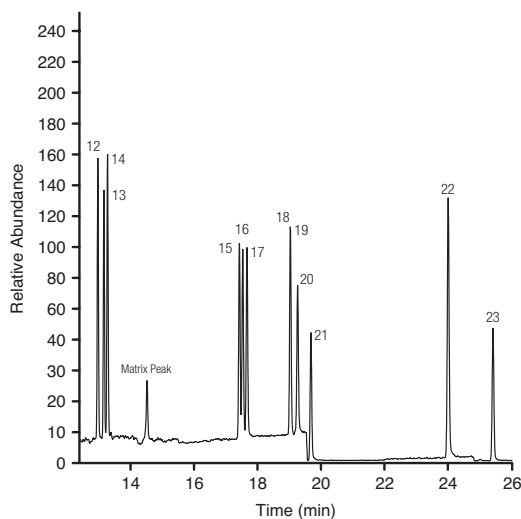


Figure 2c: Expanded SIM chromatogram of 18 PAHs in soil matrix solution at 500 ng/mL

| Peak | Compound                            | t <sub>r</sub> (min) | Linearity (R <sup>2</sup> ) | % Recovery (n=6) | %RSD (n=6) |
|------|-------------------------------------|----------------------|-----------------------------|------------------|------------|
| 1    | Napthalene-d <sub>8</sub> (IS 1)    | 4.06                 | -                           | -                | -          |
| 2    | Napthalene                          | 4.08                 | 0.9974                      | 85.0             | 0.6        |
| 3    | Acenaphthylene                      | 5.56                 | 0.9988                      | 98.2             | 1.1        |
| 4    | Acenaphthene-d <sub>10</sub> (IS 2) | 5.63                 | -                           | -                | -          |
| 5    | Acenaphthene                        | 5.66                 | 0.9992                      | 96.3             | 0.8        |
| 6    | Fluorene                            | 6.09                 | 0.9991                      | 100.2            | 1.1        |
| 7    | Phenanthrene-d <sub>10</sub> (IS 3) | 7.14                 | -                           | -                | -          |
| 8    | Phenanthrene                        | 7.15                 | 0.9992                      | 98.6             | 0.5        |
| 9    | Anthracene                          | 7.18                 | 0.9992                      | 96.7             | 0.3        |
| 10   | Fluoranthene                        | 8.98                 | 0.9985                      | 105.6            | 1.2        |
| 11   | Pyrene                              | 9.58                 | 0.9986                      | 106.7            | 1.0        |
| 12   | Benzo[a]anthracene                  | 12.98                | 0.9990                      | 99.3             | 0.4        |
| 13   | Chrysene-d <sub>12</sub> (IS 4)     | 13.16                | -                           | -                | -          |
| 14   | Chrysene                            | 13.27                | 0.9993                      | 96.9             | 0.5        |
| 15   | Benzo[b]fluoranthene                | 17.42                | 0.9993                      | 93.1             | 0.6        |
| 16   | Benzo[k]fluoranthene                | 17.54                | 0.9990                      | 93.1             | 0.6        |
| 17   | Benzo[j]fluoranthene                | 17.70                | 0.9993                      | 97.4             | 0.4        |
| 18   | Benzo[a]pyrene                      | 19.03                | 0.9990                      | 88.7             | 0.4        |
| 19   | Benzo[e]pyrene                      | 19.25                | 0.9994                      | 87.1             | 0.8        |
| 20   | Perylene-d <sub>12</sub> (IS 5)     | 19.77                | -                           | -                | -          |
| 21   | Indeno[1,2,3-cd]pyrene              | 22.03                | 0.9993                      | 91.7             | 2.2        |
| 22   | Dibenzo[a,h]anthracene              | 24.00                | 0.9993                      | 95.8             | 2.8        |
| 23   | Benzo[g,h,i]perylene                | 25.40                | 0.9990                      | 90.8             | 1.2        |

Table 3: Peak identification, linearity and recovery data for 18 PAHs in soil spiked at 1 mg/kg

## Conclusion

The QuEChERS sample preparation method provided a fast and simple approach for extracting and analyzing 18 PAHs in soil achieving high recoveries and excellent reproducibility. The QuEChERS – GC-MS method was found to be linear in the concentration range of 50 to 2000 ng/g. The TraceGOLD TG-17SilMS GC column provided good chromatographic separation for all analytes studied.

## References

1. M. Anastassiades, S.J. Lehotay, D. Stajnbaher and F.J. Schenck, J AOAC Int 86 (2003) 412.

[thermoscientific.com/chromatography](http://thermoscientific.com/chromatography)

© 2013 Thermo Fisher Scientific Inc. All rights reserved. Vespel is a registered trademark of E.I. du Pont de Nemours and Co. All other trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. This information is presented as an example of the capabilities of Thermo Fisher Scientific Inc. products. It is not intended to encourage use of these products in any manners that might infringe the intellectual property rights of others. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

**USA and Canada** +1 800 332 3331  
**France** +33 (0)1 60 92 48 34  
**Germany** +49 (0) 2423 9431 20 or 21  
**United Kingdom** +44 (0)1928 534110  
**Japan** +81 3 5826 1615

**China** +86 21 68654588 +86 10 84193588  
+86 20 83145199 800 810 5118  
**India** +91 22 6742 9494 +91 27 1766 2352  
**Australia** 1 300 735 292 (free call domestic)  
**New Zealand** 0800 933 966 (free call domestic)  
**All Other Enquiries** +44 (0) 1928 534 050

**Technical Support**  
**North America** +1 800 332 3331  
**Outside North America** +44 (0) 1928 534 440

**Thermo**  
**SCIENTIFIC**

Part of Thermo Fisher Scientific