

Whitepaper

A high-quality Raman microscope can be recognized by these 5 criteria

Damon Strom, Sonja Breuninger, Harald Fischer, Andrea Richter, Ute Schmidt, Olaf Hollricher
WITec GmbH, Lise-Meitner-Str. 6, 89081 Ulm, Germany

90 years ago Chandrasekhara Venkata Raman and Kariamanickam Srinivasa Krishnan first documented "A New Type of Secondary Radiation," which then became known as the Raman Effect [1, 2]. Raman spectroscopy is based on this effect and is used for qualitative and quantitative analysis of the chemical components and molecules of a sample. It is a nondestructive method that requires little, if any, sample preparation.

Nevertheless, Raman spectroscopy long remained a technique that was only performed in special laboratories. In recent years, however, it has been increasingly losing its outsider status. The reason for this is the development of the confocal Raman microscope, with which not only individual Raman spectra, but also complete images generated from thousands of spectra can be acquired. Through continuous development, commercially-available Raman microscopes are also becoming more user-friendly. For example, modern software interfaces guide the user through the Raman measurement and the subsequent data analysis.

There are several key factors that can be used as criteria for determining the quality of confocal Raman microscopes.

1) Speed

While in the past exposure times of minutes to hours were common for acquiring single Raman spectra, today these times are generally fractions of a second to less than one millisecond. In one second more than 1000 Raman spectra can be recorded. Thus a Raman image can be generated within a few minutes. To achieve this acquisition speed, the Raman imaging system should be equipped with optimized optics and an EMCCD camera.

High acquisition speeds are particularly important for measurements on sensitive and valuable samples in which the excitation energy must be as low as possible. Time-resolved investigations of dynamic processes can also benefit from rapid Raman spectral acquisition. Operating costs can also be reduced by shorter analysis times concurrent with increased data rates. Having a high system speed is also advantageous for time-critical work.

2) Sensitivity

The signal sensitivity of a system is critical for the quality of the results and is especially important when weak Raman signals are to be detected.

In order to achieve the best possible sensitivity, a confocal beam path, i.e. using a diaphragm aperture, must be employed to eliminate light from outside the focal plane to increase the signal-to-noise ratio. The entire Raman imaging system should also be optimized for high light throughput. This includes a spectrometer that ensures throughput of over 70 percent and is designed for measurements with low light and signal intensity. CCDs optimized for spectroscopy, which exhibit more than 90% quantum efficiency in the visible range, are most commonly used as detectors. Finally, the use of almost lossless photonic fibers ensures efficient light and signal transmission.

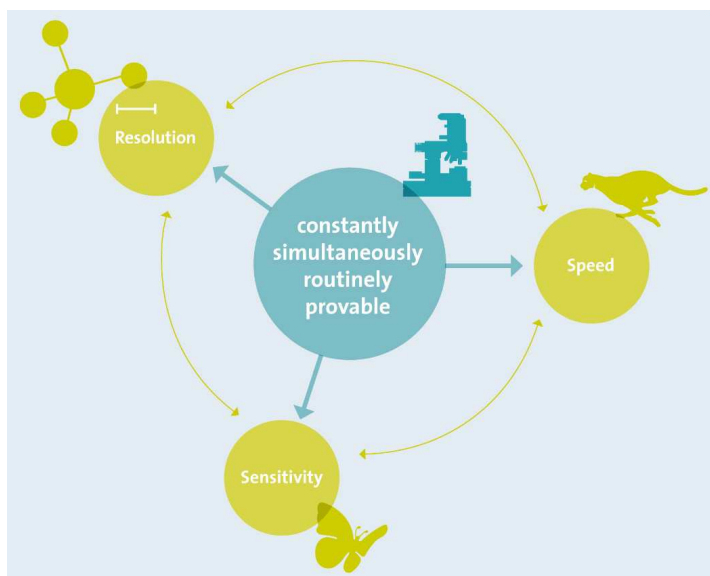


Fig. 1: Speed, sensitivity and resolution are some of the characteristics that can be used to identify a high-quality Raman microscope. These three characteristics should not be mutually exclusive. Ideally, the Raman imaging system should be configured in such a way that high-resolution images with a high signal-to-noise ratio can be acquired in a short period of time.

3) Resolution

The resolution of a Raman system is comprised of both spatial and spectral resolution. The spatial resolution includes the lateral resolution (x- and y-directions) and the depth resolution (z-direction). It is determined by the numerical aperture of the objective used and the excitation wavelength. In addition, a confocal microscope produces images with a higher contrast because the background signal is reduced. The smaller the aperture of a confocal microscope, the higher its resolution is. In a confocal Raman microscope, the lateral resolution is about 200 - 300 nm and the depth resolution below 1 μm . A confocal microscope can also create optical sections from different focal planes, which can be used with transparent samples for depth profiles and 3D images.

Spectral resolution defines the ability of a spectroscopic system to separate Raman lines near one another. Symmetric peaks in the spectrum are ensured by a spectrometer design that operates free of coma and astigmatism. The grating used, the focal length of the spectrometer, the pixel size of the CCD camera and the size of the aperture also affect the spectral resolution.

At room temperature, the width of the Raman lines is typically greater than 3 cm^{-1} , but some applications (gases, low temperature or stress analysis) may require significantly higher resolution (Fig. 1).

4) Modularity and Upgradeability

The introduction of Raman microscopy into laboratories places new demands on commercially-available systems. These requirements can sometimes appear contradictory: easy operation with diverse functionality, a wide range of applications with optimized sensitivity, low cost and high performance. In order to offer users a Raman system tailored to their individual requirements, it is particularly important that systems have a modular design that can be adapted to new conditions through being reconfigured or upgraded. A system can be optimized for specific requirements by individually combining suitable lasers, filters, lenses, spectrometers and detectors. With such a customized Raman imaging system the user is able to obtain meaningful Raman images, perform 3D volume scans and create depth profiles.

5) Combinability

Confocal Raman microscopy can be combined with other microscopy techniques. By using different methods and correlating the data, the user attains a more comprehensive understanding of the sample. Common examples of correlative microscopy techniques are Raman-AFM, Raman-SNOM and Raman-SEM. In order to correlate the data of these disparate technologies, the exact same sample location must be examined by each approach. If different instruments are to be used, finding this sample location can be very difficult and time-consuming. This is made much easier with a hybrid system that combines the different analysis methods in one instrument so that the sample can remain in place during all measurements.

Some applications of correlative Raman microscopy are presented below.

Raman and Profilometry

For Raman microscopy, most samples do not need to be treated, stained or otherwise prepared prior to measurement. The combination of a confocal Raman microscope with a profilometer module for focus stabilization allows rough or inclined surfaces to be examined [3, 4]. During Raman analysis, the examination area is kept constantly in focus by the simultaneously-acquired profilometry data. This also compensates for thermal shifts and enables long-term measurements. The application example in Fig. 2 shows the analysis of a microstructured silicon sample. The chemical image of the Raman measurement was overlaid onto the topographic profile measurement.

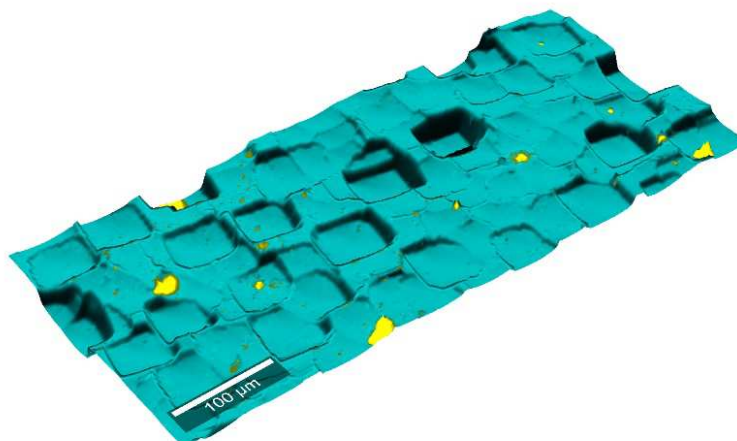


Fig. 2: Topographic Raman image of a silicon microstructure.

Raman and Fluorescence

Fluorescence microscopy has been a widespread imaging method for the analysis of biological cells and organisms for decades. Samples are stained with fluorescent dyes or organisms are genetically engineered to express fluorescent proteins. The fluorescence signal is usually much stronger than the Raman signal. Nevertheless, correlative Raman fluorescence measurements are possible with an appropriate system. Fig. 3 shows a Raman fluorescence image of a live cell culture of eukaryotic cells. An inverted confocal Raman microscope was used to examine the cells in their aqueous cell culture medium in the Petri dish. The cell nuclei were stained with the fluorescent dye DAPI. An excitation wavelength of 532 nm was used for the Raman measurement. An image with $50 \times 40 \mu\text{m}^2$ and 150×120 pixels was acquired. A Raman spectrum was recorded at each pixel. The recording time was 0.2 s/spectrum. In the correlative Raman fluorescence image, the nuclei are shown in blue (recorded with fluorescence microscopy), the nucleoli in green and the endoplasmic reticula in red (recorded with Raman microscopy). The corresponding Raman spectra are shown in the same colors.

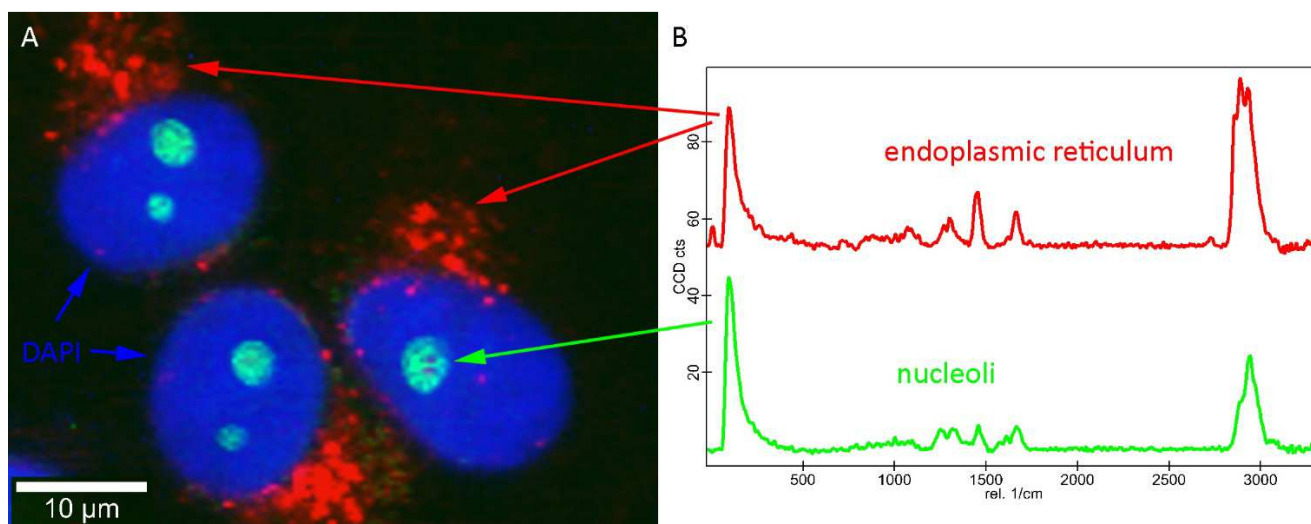


Fig. 3: A: Correlative Raman fluorescence image of primate cells in cell culture. Blue: Nuclei recorded by fluorescence microscopy; Red: Endoplasmic reticula, Green: Nucleoli recorded by Raman microscopy. B: Associated Raman spectra.

Raman and AFM

The combination of Raman microscopy, which provides information about the type and distribution of molecules in a sample, and the high-resolution AFM technique, which determines the surface characteristics of a sample, enables the visualization of both chemical and morphological properties. Here the analysis of a 1:1:1 mixture of polystyrene (PS), 2-ethylhexyl acrylate (EHA) and styrene-butadiene rubber (SBR) is shown. For this, a correlative Raman-AFM microscope was used, in which Raman microscopy and AFM technologies are fully integrated.

The measurement with AFM in intermittent AC mode documents the topography of the polymer mixture (Fig. 4A). The simultaneously-recorded phase image (Fig. 4B) provides information on the viscosity and

elasticity of the individual components of the polymer mixture. The confocal Raman image (Fig. 4C) shows that PS (red) and EHA (green) are present separately. SBR (purple) partly mixes with EHA (mixture shown in blue). By correlating the Raman image with the AFM image, the chemical information can be linked to the structural information (Fig. 4D).

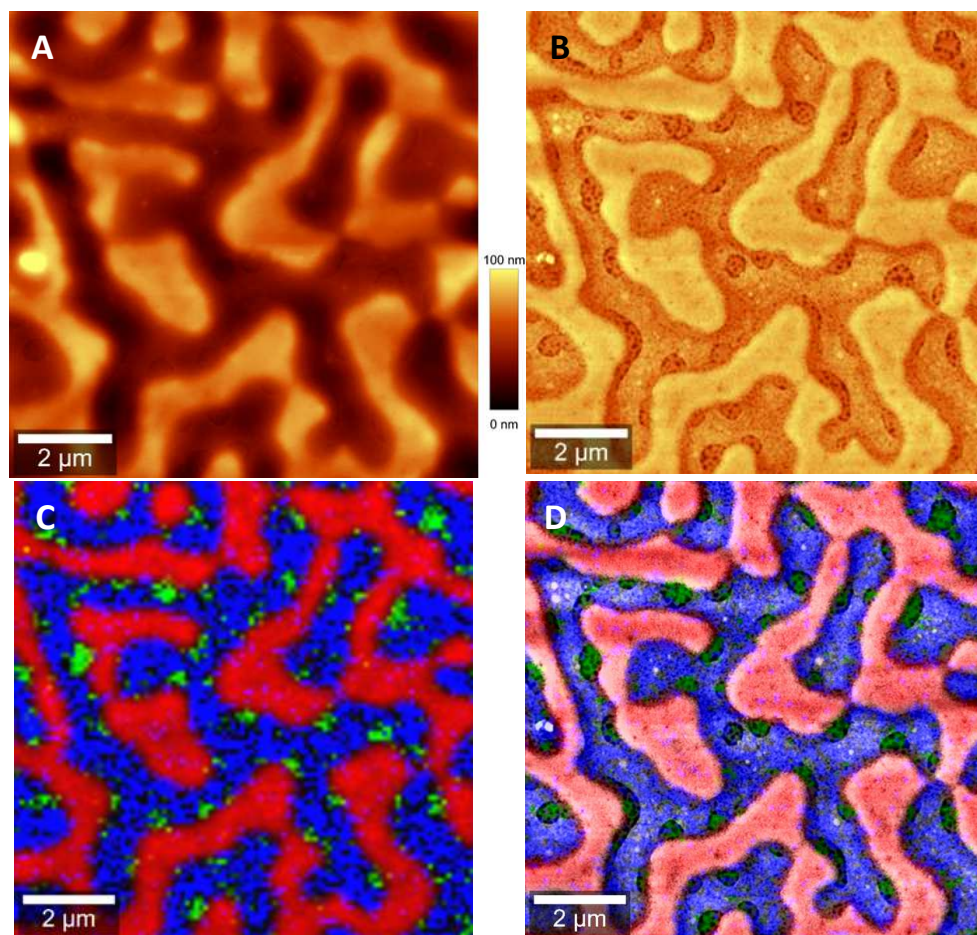


Fig. 4 : Correlative, high-resolution AFM-Raman microscopy of a 1:1:1 mixture of polystyrene (PS), 2-ethylhexyl acrylate (EHA) and styrene-butadiene rubber (SBR). Both measurement technologies are combined in the WITec alpha300 RA microscope. A: The topography of the polymer mixture was determined with AFM in AC mode. B: The phase of the AFM image shows the fine structure of the compound. C: A color-coded, confocal Raman image, generated from the Raman spectra, shows the distribution of the polymers. PS (red), EHA (green). SBR (purple), SBR-EHA mixture (blue). D: In the correlative Raman-AFM image the topography and distribution of the different polymers can be visualized.

Raman and SEM

Scanning electron microscopy (SEM) is a well-established method for structural surface analysis. By combining Raman imaging with SEM in a correlative microscope, it is possible to combine results of SEM structural analysis with chemical and molecular information from confocal Raman microscopy [5]. The sample is placed in the vacuum chamber of the electron microscope. Both analysis methods are then carried out automatically at the same sample location. The obtained SEM and Raman images can then be superimposed.

In Fig. 5 a structure several atoms in thickness comprised of graphene layers was analyzed by correlative Raman-SEM microscopy. The Raman image consists of 22,500 spectra with 50 ms recording time per spectrum. While in the SEM image the contrast between the substrate and the graphene flake is visible, in the Raman image the number of graphene layers and their different orientations can be analyzed. This is not possible with SEM alone.

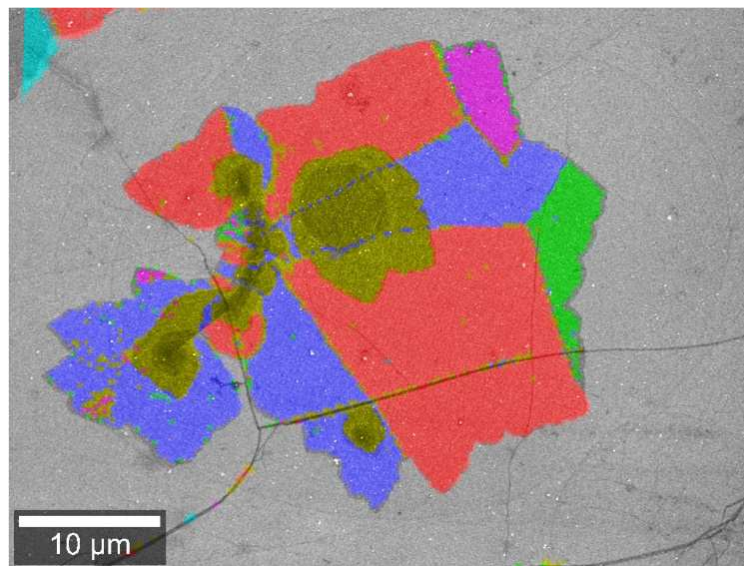


Fig. 5: Correlative RISE (Raman-SEM) microscopy image of a multilayer graphene flake. The different colors show folds and orientations in the graphene and could be identified by Raman analysis.

Raman Particle Identification and Characterization

High-resolution investigations of particles are of great interest in many fields of application such as environmental science, pharmaceutical research and many others. Combining a particle analysis tool with the fast, label-free and nondestructive Raman imaging technique makes it possible to find, classify, and identify particles in a sample according to their size, shape and chemical characteristics. The physical and molecular attributes of the particles in a sample may be correlated and qualitatively and quantitatively evaluated. Fig. 6 shows the results of particle analysis carried out on a cosmetic peeling cream sample. Fig. 6A shows the overlay of an optical bright field microscope image with the corresponding confocal Raman image. Particles are identified according to their size and shape and further characterized by their molecular properties through confocal Raman imaging. The chemical analysis revealed anatase and boron nitride particles in an oil matrix (Raman spectra shown in Fig. 6B). Further evaluation of the results determines the quantitative prevalence of the molecular sample components in the particles (Fig. 6C) and also the distribution of chemical compounds correlated to particle size (Fig. 6D). In extended analyses the chemical characteristics of particles could also be linked to parameters such as area, perimeter, bounding box, Feret diameter, aspect ratio, equivalent diameter and many others. This illustrates the potential for comprehensive investigations of particles in many fields of application.

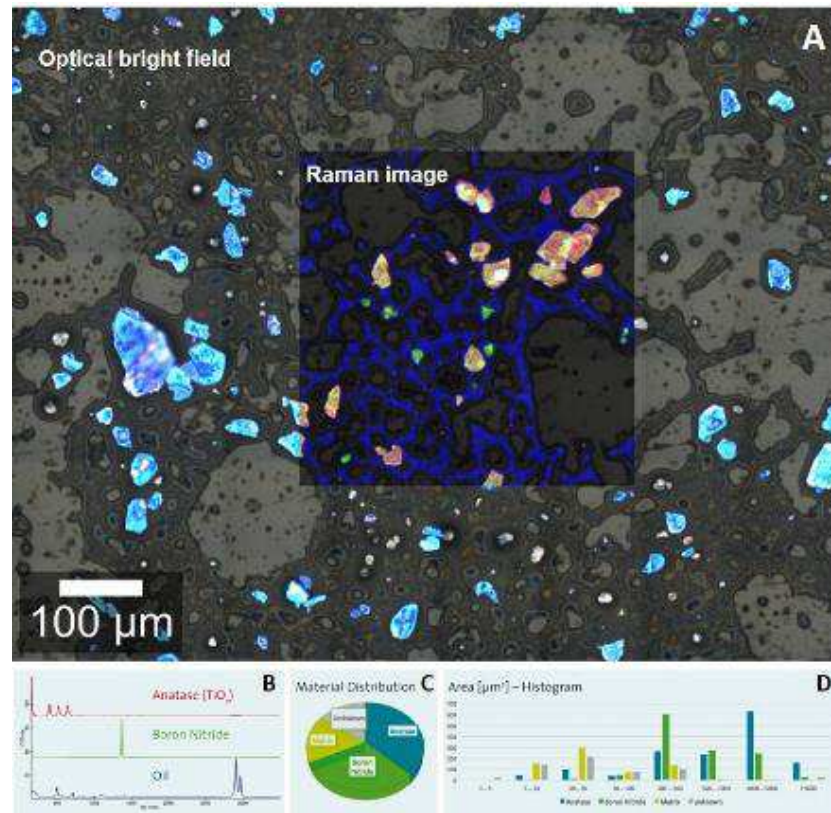


Fig. 6: Particles in a cosmetic peeling cream sample. A: Optical bright field image overlaid with the confocal Raman image. B: Corresponding Raman spectra of the molecular components in the sample. C: Pie chart of the quantitative compound distribution in the sample. D: Graphical representation of the correlation between chemical characteristics and particle size.

Literature:

- [1] C.V. Raman, K.S. Krishnan, *Nature* 121, 501 (1928)
- [2] J. Toporski, T. Dieing and O. Hollricher (eds.), *Confocal Raman Microscopy*, 2nd Edition, Springer Series in Surface Sciences (66), Springer International Publishing (2018)
- [3] A. Masic, J. C. Weaver, *Journal of Structural Biology* 189, 269-275 (2015)
- [4] B. Kann, M. Windbergs, *The AAPS Journal* 15, 505-510 (2013)
- [5] G. Wille, C. Lerouge, U. Schmidt, *Journal of Microscopy* 270, 309-317 (2018)



Contact:

Harald Fischer
Marketing Director
Harald.Fischer@witec.de

WITec GmbH
Lise-Meitner-Str. 6
89081 Ulm
Germany

Tel.: +49 (0) 731 140 70-0
Fax: +49 (0) 731 140 70-200
www.witec.de
info@witec.de

About WITec:

WITec GmbH pioneered 3D Raman imaging and correlative microscopy and continues to lead the industry with a unique product portfolio that offers speed, sensitivity and resolution without compromise. Raman, AFM, SNOM and SEM (RISE) microscopes and select combinations thereof can be configured for specific challenges in chemical and structural characterization. WITec's headquarters is located in Ulm, Germany, where all WITec products are developed and produced. Branch offices in USA, Japan, Singapore, China, France and Spain provide a worldwide sales and support network. WITec's enduring success and innovative product range have been recognized by a series of significant industry awards. Further information can be found at www.witec.de